

輝光放電機 *Glow Discharge System*

PELCO easiGlow 91000 注意事項 PELCO easiGlow 91000 Precautions

1.



打開電源
Turn on Power

2.



向上輕拿玻璃蓋、勿撞中間金屬！
CAREFULLY move the glass chamber
upward. Don't hit the metal part.



輕放到旁邊桌上
Gently put it on the
table aside.

輝光放電機 *Glow Discharge System*

3.

將網格放上載玻片，並放入樣品凹槽中。

Place grids on a slide glass and put it onto the groove.



4.



小心蓋回上蓋，紅色墊片須對準下方黑色密封條。

Carefully put the chamber back and check the red gasket alignment with the black O-ring seal.

5. 時間預設為 20 秒，請觸控螢幕上 **AUTO RUN** 的開始輝光放電。如需輝光放電 20 秒以上，請重複此步驟。

Default time is 20 seconds, hit **AUTO RUN** on the touch screen to start glow discharge. If need more than 20 seconds, please repeat this step.

Default Setting:

- SET: 20mA
- GLOW: 20s



輝光放電機 *Glow Discharge System*



6. 運轉結束時，須等洩壓完成後再小心開啟上蓋取出網格。

After finishing the glow discharge **and venting**, open the chamber carefully and take the grids out.

7. 放回上蓋，再直接關閉電源。

Replace the chamber back, then directly **Turn Off** the power.