



CEM3 (Talos Arctica) Operational Safety Precaution 操作安全注意事項！

TEM User Interface (Monitor 1)

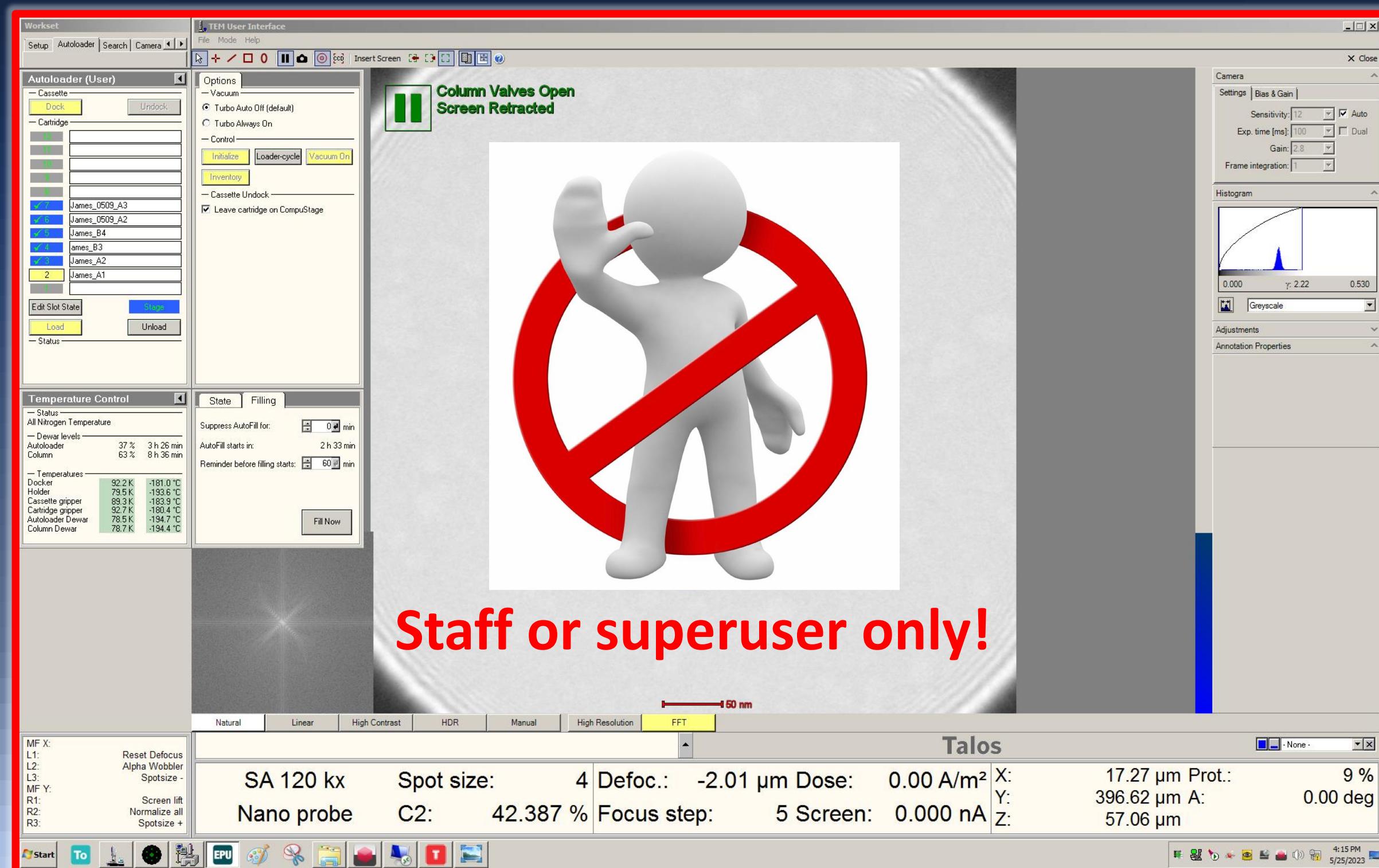

EPU (Monitor 2)

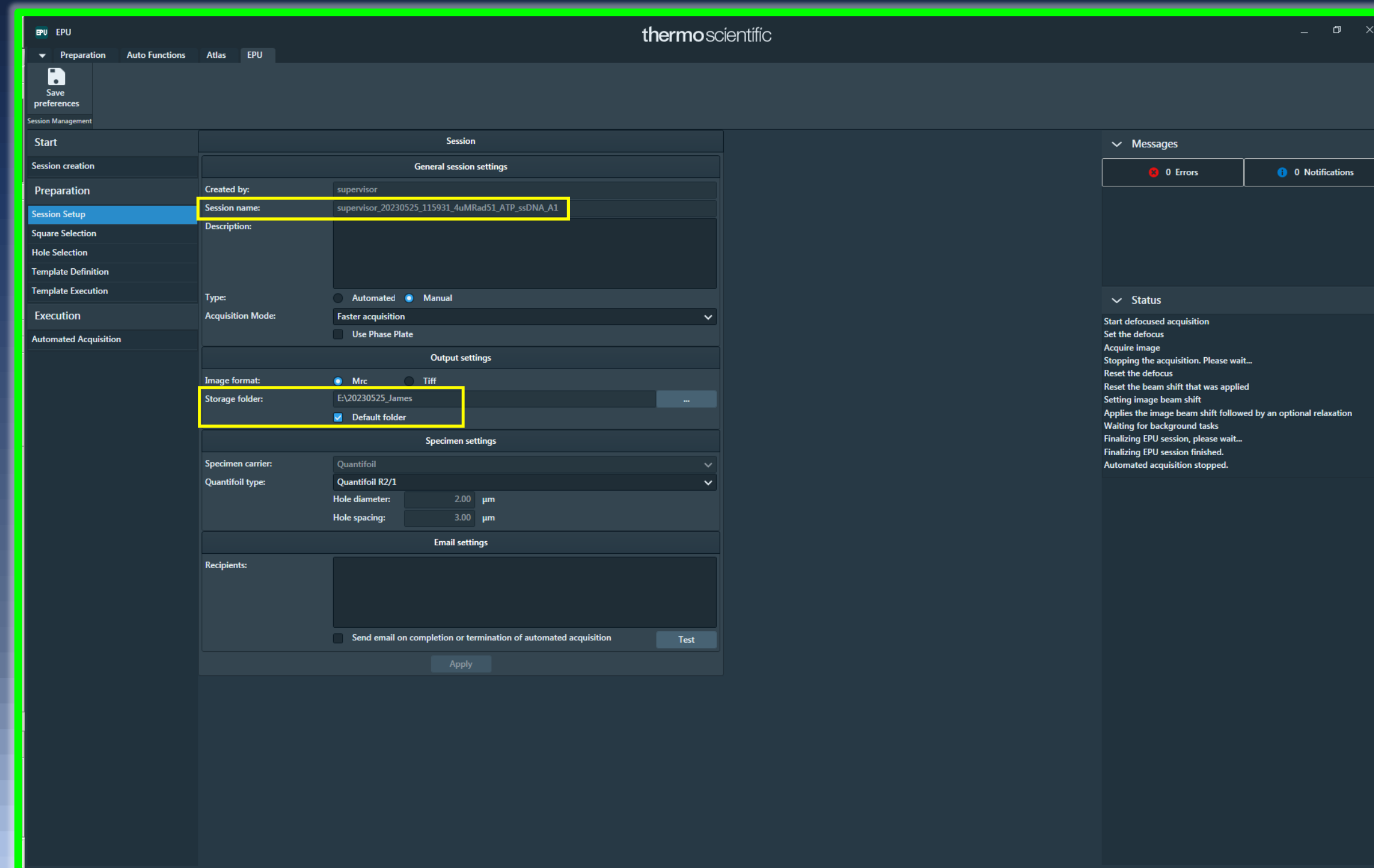

Session Setup

- Session name: **Supervisor_xxx_sample_name**
Sample Naming (樣品命名)
ONLY letters, numbers and underline are allowed for sample naming.
The specific symbol can cause the PC to CRASH!
If a CRASH happened, instrument usage fees will still be charged!
樣品命名只准許輸入英文字母、數字及底線。
使用特殊符號會造成當機！當機時間仍計入收費時段，上機費用照收！
- Storage folder: **E:\date_name**
Please select **E** drive to create your own file folder.
檔案只能存至 **E** drive。

Overall CEM3 operation is using **EPU**, **NOT need to use the TEM User Interface**.

Contact CEM3 staff if you have any problems, thank you!

CEM3操作全程都使用EPU頁面，毋需用到TEM User Interface頁面，任何問題請洽工作人員，謝謝！


If you fail to follow the CEM3 operating procedures and rules, the instrument usage fees will still be charged !

操作不當，上機費用照收！


Simple Trouble Shooting:

For atlas screening, if you see the error message below:

掃Atlas時，如果遇到以下錯誤訊息：

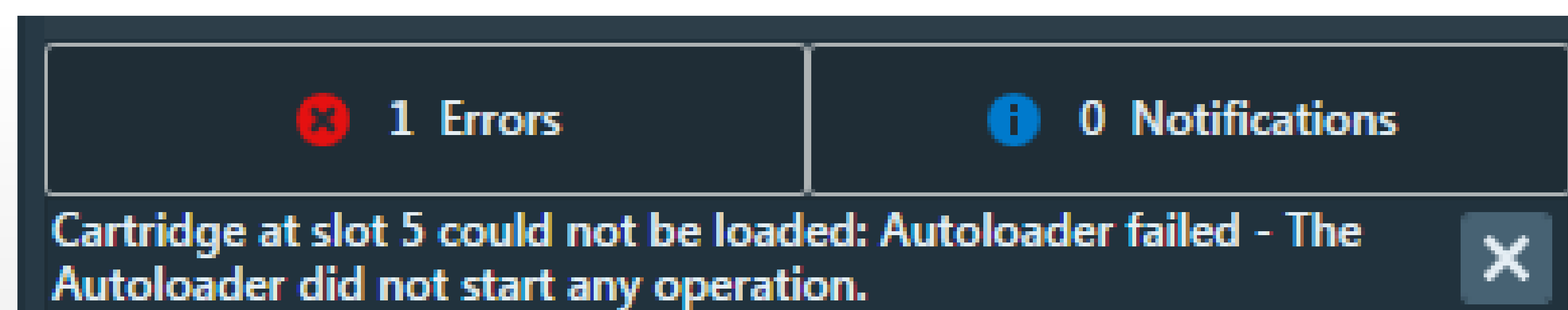

Please wait for 2 mins and click **Reset Selected**, then click **Start** again.
請等待2分鐘後按 **Reset Selected**，再按 **Start** 重新掃Atlas。

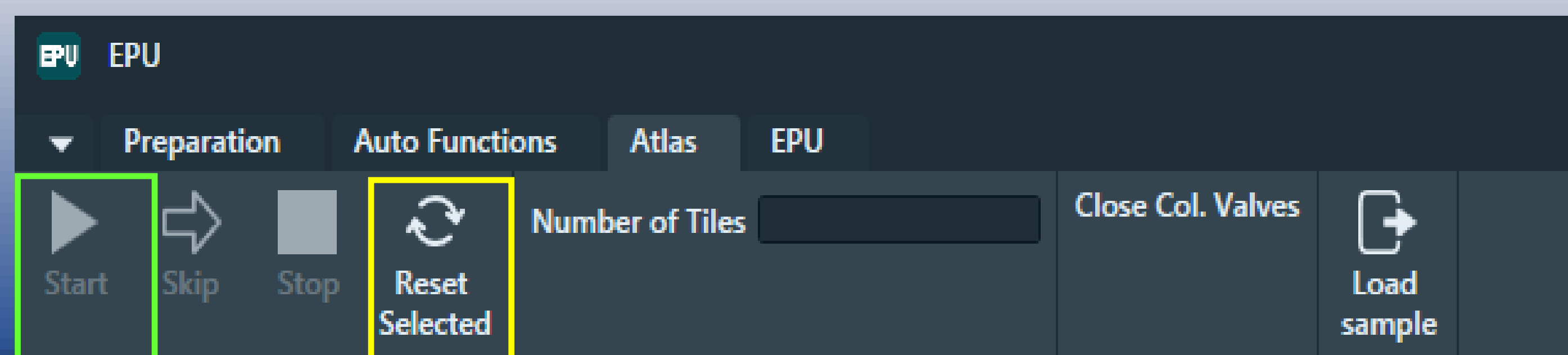

CEM3 Time Slots:

1. Morning session : 10:00 – 14:00
2. Afternoon session : 14:00 – 18:00

For afternoon session, if you want to collect data overnight, please finish the hole selection setup before 17:30 and inform CEM3 staff. 下午時段之使用者，若需要收整晚，請在 17:30 前選好洞後告知CEM3儀管人員。

Other Reminders:

1. Deliver your grids to ASCEM at least **2 business days before machine time** and submit your **grids storage information form** to CEM3 staff.
上機兩個工作天前送樣，遞交樣品放置單。
2. Record machine usage time in **the CEM3 logbook** before you leave. 記錄上機時間。
3. To recycle the screened grids, please write a note in both your **grids storage information form** and **the CEM3 logbook**. 網格回收，須同時註記在樣品放置單及記錄本。
4. To cancel the CEM3 reservation, please contact CEM3 manager at least **48 hours in advance**. 預約取消，最晚須於48小時前通知 CEM3儀管人員。

CEM 3 Staff :

1. Dr. Irene Deli 李宥慧 博士 (email : deli24@gate.sinica.edu.tw)
2. Ms. Yu-Hua Chang 張與華 小姐