




ASCEM Vitrobot 3 Upgrade Reminders

Ver. 25.2.2025 Kondo

1


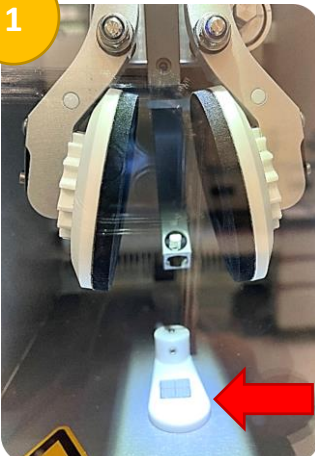


新版螢幕請直接用手指觸控操作！

In the new version,
the screen can be operated directly via touch input!

開機程序完成後，舊版的檔板狀態為開啟，新版則維持關閉狀態。
After the startup process is completed, the shutter in the original version remained open, whereas in the new version, it stays closed.

2

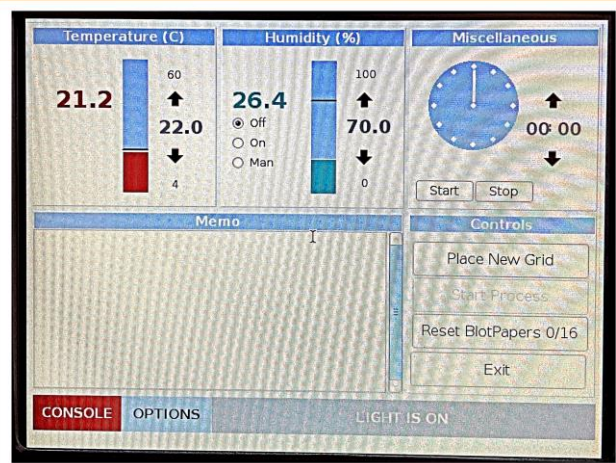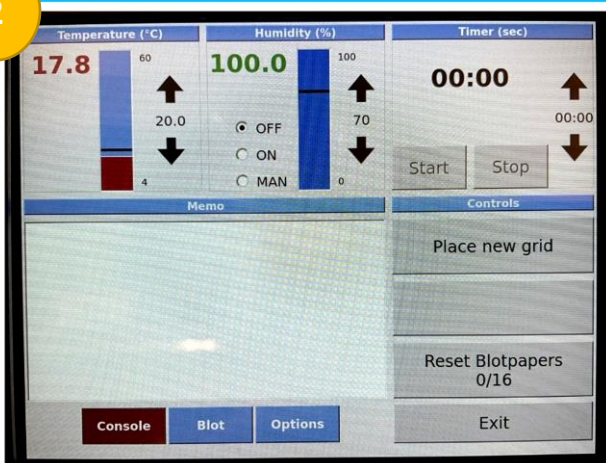



以下依序說明 **Console**、**Blot**、**Options** 的新舊版差異，新版操作介面皆如左圖所示。

- **Console**：右上角新增 Timer 功能，其餘操作與舊版一致。

Below are the key differences between the old and new versions for **Console**, **Blot**, and **Options**.
The new operation interface appears as shown in the left images.

- **Console**: A Timer function has been added in the top right corner, while all other operations remain unchanged from the previous version.

3

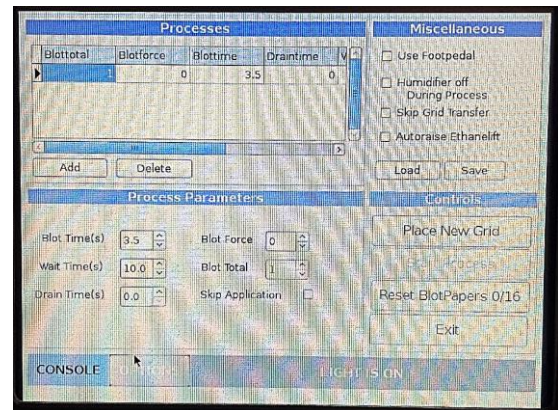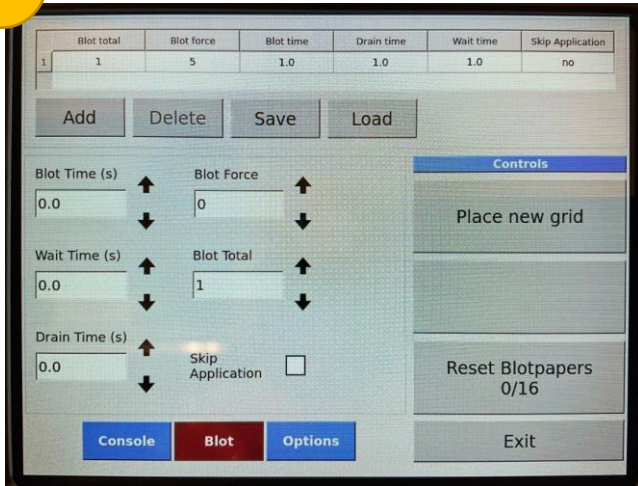



- **Blot:** 參數設定方式與舊版一致。參數設定完成後，請務必點擊 **Add** 儲存，執行 **Blot** 時將按照最上方表格中的參數執行（注意：如果參數設定未點擊 **Add**，系統將以 0 秒執行）。

新功能 若表格中輸入兩組以上參數，則可執行 **Multiple Blotting**。

- **Blot:** The parameter settings remain unchanged from the previous version. Please ensure to click **Add** after configuring the parameters. When executing **Blot**, the system will follow the parameters in the topmost table. (Note: If **Add** is not clicked, the system will default to 0 second.)

New Feature If more than two sets of parameters are entered in the table, **Multiple Blotting** can be performed.

4

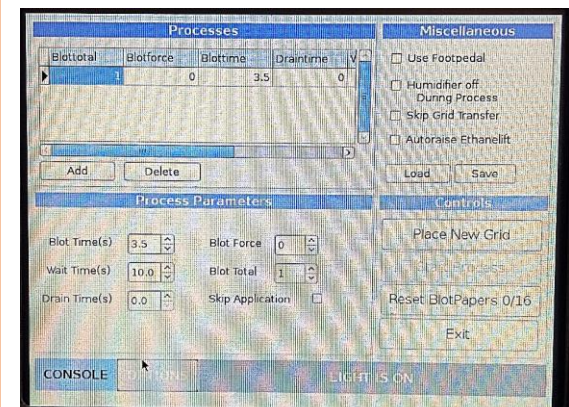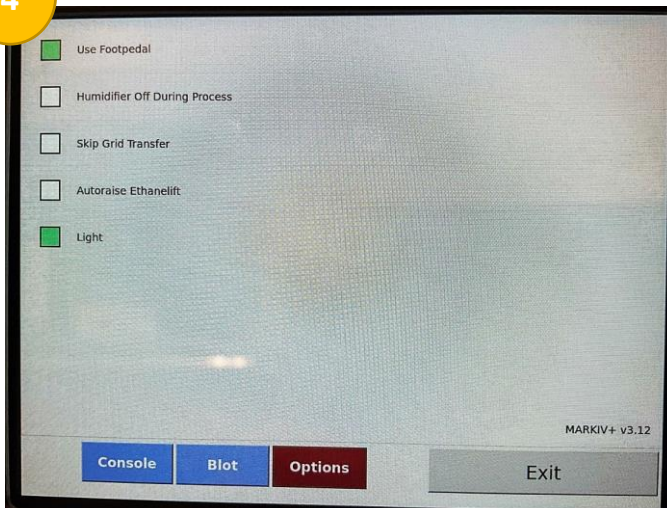

- **Options:** 參數設定方式與舊版一致。此畫面在舊版中顯示於 **Options** 介面的右上角。
- **Options:** The parameter settings remain unchanged from the previous version. In the old version, this part is located in the top right corner of the **Options** page.